

Nitrogen/Protein Determination in Animal feed by the Thermo Scientific FLASH 2000 Analyzer using Argon as Carrier Gas

Dr. Liliana Krotz, OEA Product Specialist, Walter Galotta, OEA Product Support Engineer, and Dr. Guido Giazzi, OEA Product Manager, Thermo Fisher Scientific, Milan, Italy

Key Words

Animal feed, argon, combustion, nitrogen, protein

Goal

This application note presents data on nitrogen/protein determination in reference materials, raw and final animal products with different nitrogen concentrations to show the performance of the system using argon as an alternative carrier gas and show the reproducibility of the results obtained.

One of the most important nutrients in animal nutrition is protein; the intake of which provides the body with a resource for energy, growth and production of muscles, hormones, milk, wool etc.

Protein is a compound of amino acids. Essential amino acids are those which the body cannot create and must therefore be taken in as food. Essential amino acids differ between species, and as such the level of proteins/amino acids in animal food must exist in specific and correct levels to achieve the nutritional quality of the finished product. They must provide a balanced level of protein to prevent excess or avoid a deficiency in amino acids. This can be monitored through the precise and accurate determination of the levels of nitrogen in the product.



Figure 1. FLASH 2000 Elemental Analyzer



For this reason, modern advances in instrumentation for the determination of nitrogen have greatly improved the capabilities of the combustion method, making it faster, safer and more reliable than the traditional Kjeldahl method. The Dumas (combustion) method has been approved and adopted by the Association of Official Analytical Chemists (AOAC Official Method 990.03. Protein crude in Animal Feed 4.2.08).

The supply of helium over the last few years has been beset with problems, namely world-wide shortages and subsequent price increases. It has therefore been necessary to test an alternative gas, argon, which is readily available.

The Thermo Scientific™ FLASH 2000 analyzer, based on the dynamic flash combustion of the sample, copes effortlessly with a wide array of laboratory requirements, such as accuracy, day by day reproducibility and high sample throughput.

Thermo
SCIENTIFIC

Methods

Samples are weighed in tin capsules and introduced into the combustion reactor via the Thermo Scientific™ MAS™ 200R autosampler together with the proper amount of oxygen. After combustion, the resultant gases are carried by an argon flow to a second reactor filled with copper, then swept through CO₂ and H₂O traps, a GC column and finally detected by a thermal conductivity detector (TCD). The analytical configuration as well as the TCD detector are the same as those used with helium as the carrier gas (see Figure 2).

A complete report is automatically generated by the Thermo Scientific™ Eager Xperience data handling software and displayed at the end of the analysis. The Eager Xperience software provides a new option AGO (Argon Gas Option) through which modifies the argon carrier flow during the run to optimize the analysis.

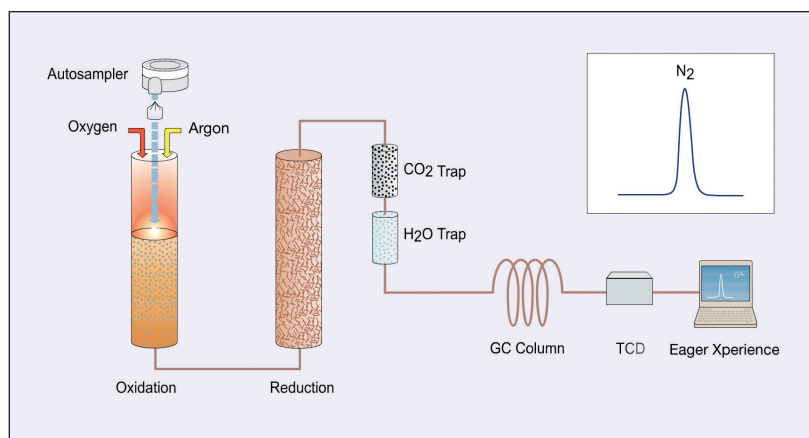


Figure 2 – FLASH 2000 Nitrogen configuration

Analytical conditions

Combustion Furnace Temperature:	950 °C
Reduction Furnace Temperature:	840 °C
Oven Temperature:	50 °C (GC column inside the oven)
Argon Carrier Flow:	60 ml/min
Argon reference Flow:	60 ml/min
Oxygen Flow:	300 ml/min
Sample Delay:	10 sec
Run time:	10 mins

Note: The oxygen amount necessary for the complete combustion of samples is calculated automatically by the OxyTune function present in the Eager Xperience software.

Official methods requirements for Protein determination by combustion

AOAC (Method 990.03) indicates that a suitable fineness of grind must be determined (for each different material analyzed) to achieve precision that gives RSD of $\leq 2\%$ for 10 successive determinations of nitrogen.

Different reference materials, raw and final animal feed products were chosen to demonstrate the availability of the method for all range of protein content without matrix effect.

Tables 1 and 2 show the sample information and the reproducibility of 10 consecutive analyses of BIPEA Reference Materials using a sample weight of approximately 100 – 120 mg for feed for sow and hyperproteic powder, and 100 – 140 mg for alfalfa. The materials were characterized through a laboratory intercomparison using Kjeldahl and Combustion methods. The protein factor used to calculate the protein content was the default value 6.25 present in the Eager Xperience software. Instrument calibration was performed with Aspartic acid (10.52 %N) standard using K factor as calibration method. The data obtained are inside the tolerance declared in the relative certificates.

Table 1 – Bipea sample information available

Sample	Moisture	Fat	Carbohydrate	Kjeldahl Protein		Combustion Protein	
	%	%	%	Av. %	Tolerance	Av. %	Tolerance
Bipea - Feed for Sow 3/2009	9.8	2.8	48.7	16.0	0.6	16.2	0.6
Bipea - Dehydrated Alfalfa 3/2009	7.7	---	29.3	14.8	0.6	15.1	0.6
Bipea - Hyperproteic Powder 1/2008	---	0.8	---	85.4	3.4	86.4	3.5

Table 2 – Reproducibility of Nitrogen/Protein determination in Bipea Reference Materials

Sample	Bipea – Feed for Sow		Bipea – Dehydrated Alfalfa		Bipea - Hyperproteic Powder	
	N %	Protein %	N %	Protein %	N %	Protein %
	2.67	16.71	2.47	15.46	13.53	84.59
	2.67	16.67	2.49	15.57	13.56	84.74
	2.66	16.63	2.45	15.34	13.67	85.42
	2.60	16.22	2.41	15.04	13.62	85.10
	2.63	16.41	2.47	15.41	13.69	85.57
	2.67	16.72	2.48	15.49	13.66	85.36
	2.67	16.66	2.37	14.81	13.60	85.02
	2.67	16.71	2.44	15.27	13.63	85.21
	2.61	16.32	2.38	14.87	13.71	85.72
	2.65	16.55	2.37	14.79	13.67	85.43
Average %	2.65	16.56	2.43	15.21	13.63	85.22
RSD %	1.02	1.09	1.93	1.97	0.42	0.42

Table 3 shows the Nitrogen/Protein data of different animal feed samples using argon and helium carrier gas including the range of weight used for analysis. The protein factor used was 6.25. No significant difference was observed.

Sample	FLASH 2000 using Argon carrier gas					FLASH 2000 using Helium carrier gas				
Name	Weight (mg)	N %	RSD %	Protein %	RSD %	Weight (mg)	N %	RSD %	Protein %	RSD %
Wheat flour	130 - 140	1.91 1.93 1.94	0.79	11.93 12.05 12.14	0.87	160 - 200	1.92 1.93 1.90	0.79	11.98 12.06 11.88	0.75
Maize 1	130 - 140	1.25 1.25 1.26	0.46	7.84 7.83 7.87	0.27	150 - 200	1.26 1.26 1.27	0.46	7.88 7.93 7.91	0.32
Maize 2	130 - 140	1.48 1.49 1.49	0.39	9.24 9.29 9.29	0.31	150 - 200	1.50 1.50 1.49	0.38	9.41 9.39 9.34	0.38
DDGS	120 - 130	4.79 4.78 4.83	0.55	29.92 29.85 30.21	0.64	150 - 160	4.79 4.80 4.77	0.31	29.92 30.03 29.80	0.38
Sunflower 1	100 - 120	6.02 6.08 6.09	0.62	37.62 38.01 38.08	0.65	170 - 200	6.10 6.07 6.09	0.25	38.15 37.91 38.07	0.31
Sunflower 2	130 - 140	5.69 5.75 5.71	0.53	35.54 35.95 35.68	0.58	170 - 200	5.65 5.61 5.64	0.37	35.31 35.05 35.22	0.37
Soya	130 - 150	8.11 8.08 8.00	0.71	50.66 50.51 49.99	0.70	170 - 180	7.98 8.03 8.01	0.31	49.87 50.21 50.06	0.34
Gluten 1	130 - 140	9.31 9.27 9.28	0.22	58.20 57.92 58.02	0.24	170 - 180	9.34 9.33 9.31	0.16	58.36 58.33 58.21	0.14
Gluten 2	130 - 140	2.91 2.98 2.96	0.38	58.18 58.44 57.98	0.40	170 - 180	9.32 9.37 9.36	0.28	58.28 58.55 58.50	0.25
Poultry feed 1	120 - 140	2.91 2.98 2.96	1.22	18.18 18.62 18.48	1.22	150 - 160	2.95 2.96 2.90	1.09	18.43 18.53 18.13	1.13
Poultry feed 2	130 - 140	4.31 4.39 4.30	1.14	26.96 27.43 26.87	1.11	150 - 160	4.37 4.35 4.42	0.82	27.30 27.17 27.65	0.91

Conclusions

Good repeatability, accuracy and precision was obtained with the FLASH 2000 nitrogen analyzer using argon as carrier gas to analyze nitrogen in a wide range from low to high content without matrix effect.

No memory effect was observed when changing the type of sample, indicating complete combustion and detection of the element. Samples used were chosen for their high content of protein.

The nitrogen/protein data obtained of the BIPEA Reference Materials are inside the tolerance declared in relative certificates and the RSD % obtained was less than 2% according to Official Methods Performance Requirements.

www.thermoscientific.com

©2014 Thermo Fisher Scientific Inc. All rights reserved. BIPEA is a trademark of the Bureau Interprofessionnel des Etudes Analytiques (International Bureau for Analytical Studies). AOAC is a trademark of The Association of Official Analytical Chemists. ISO is a trademark of the International Standards Organization. All other trademarks are the property of Thermo Fisher Scientific and its subsidiaries. This information is presented as an example of the capabilities of Thermo Fisher Scientific products. It is not intended to encourage use of these products in any manners that might infringe the intellectual property rights of others. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

COMPANY WITH QUALITY MANAGEMENT
SYSTEM CERTIFIED BY DNV
= ISO 9001:2008 =

Thermo Fisher Scientific S.p.A. Milan, Italy
is ISO 9001:2008 Certified.

Africa +43 1 333 50 34 0
Australia +61 3 9757 4300
Austria +43 810 282 206
Belgium +32 53 73 42 41
Canada +1 800 530 8447
China 800 810 5118 (free call domestic)
400 650 5118

Denmark +45 70 23 62 60
Europe-Other +43 1 333 50 34 0
Finland +358 9 3291 0200
France +33 1 60 92 48 00
Germany +49 6103 408 1014
India +91 22 6742 9494
Italy +39 02 950 591

Japan +81 45 453 9100
Latin America +1 561 688 8700
Middle East +43 1 333 50 34 0
Netherlands +31 76 579 55 55
New Zealand +64 9 980 6700
Norway +46 8 556 468 00
Russia/CIS +43 1 333 50 34 0

Singapore +65 6289 1190
Spain +34 914 845 965
Sweden +46 8 556 468 00
Switzerland +41 61 716 77 00
UK +44 1442 233555
USA +1 800 532 4752

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific